

**Specification Sheet**

**Product:      Drying Buffer/Stabilizer, biotin-free BSA based,  
                  ready-to-use**

**Code: #DBS-B**

**Protein concentration :**           10 mg/ml

**Buffer base:**                    0,15M PBS pH 7,3

**Preservative:**                    BND, 550 ppm

**Lot:**                                765123

**Storage:**                         +2°C ~ +8°C

**Expiration date:**                05/2014

**Recommended use:**                ready-to-use reagent for  
preparing dried non porous and  
porous immunosorbents that shall be stable real-time at  
room/ambient temperature

Warm up to room temperature prior to use

It is suggested that e.g. for ELISA work DBS-C will be applied AFTER washing of the ImmunoPlate with the wash buffer following antibody and/or antigen adsorption, i.e. BEFORE drying. Do not wash DBS-C out of the plate, , just aspirate it as completely as possible immediately before drying.